

## 1 Background

As the need for multiple biomarker assessment in immunology has become more clear, multiplex fluorescent immunohistochemistry (IHC) techniques have become integral to immune oncology research.

Multispectral approaches to IHC improve quantitative performance and signal specificity by ensuring removal of autofluorescence signals and bleed-through between spectrally-adjacent dyes.

We have recently demonstrated a novel high-throughput scanning approach (MOTIF™) that allows acquisition of a 7-color multispectral slide scan in 5–10 minutes (1.5x1.5 cm).

Here, we demonstrate the spatial measurements that are made possible by this high-throughput, translational workflow. These measurements span the scales related to tumor biology from the distance between nearest cell neighbors to the extent of the invasive margin.

## 2 Methods


### Phenoptics™ Workflow

Formalin-fixed paraffin-embedded (FFPE) samples of primary tumors were stained using Opal™ reagents on a Leica BOND RX. Sections from several primary tumor types were stained against 5 key markers in the field of immunology (CD8, FoxP3, PD-1, PD-L1, CD68) along with a tumor marker (PanCK, PAX8, or Sox10+S100 cocktail) and DAPI counterstain, resulting in 7 colors on each slide.

Conventional and MOTIF™ multispectral scans (0.5 μm/pixel; 2.8 – 7.4 cm²) were acquired on Vectra Polaris® (9 – 26 min scan times). Scans were viewed with a Phenochart™ version modified to divide analysis regions into larger fields to reduce batch analysis times. Cell segmentation and phenotyping were performed in inForm®. Batch results from inForm® were merged and consolidated with the phenoptReports package<sup>1</sup> in RStudio<sup>2</sup> using R 3.5.2<sup>3</sup>. Tumor margin annotations were drawn and exported with QuPath<sup>4</sup>. Heatmaps, histograms, and density plots were generated from consolidated inForm® results with R scripts under development.

### Multiplex Staining

Opal™ reagents allow multiplex IHC staining of up to 8 antigens with signal amplification and any combination of mouse and/or rabbit primary antibodies.

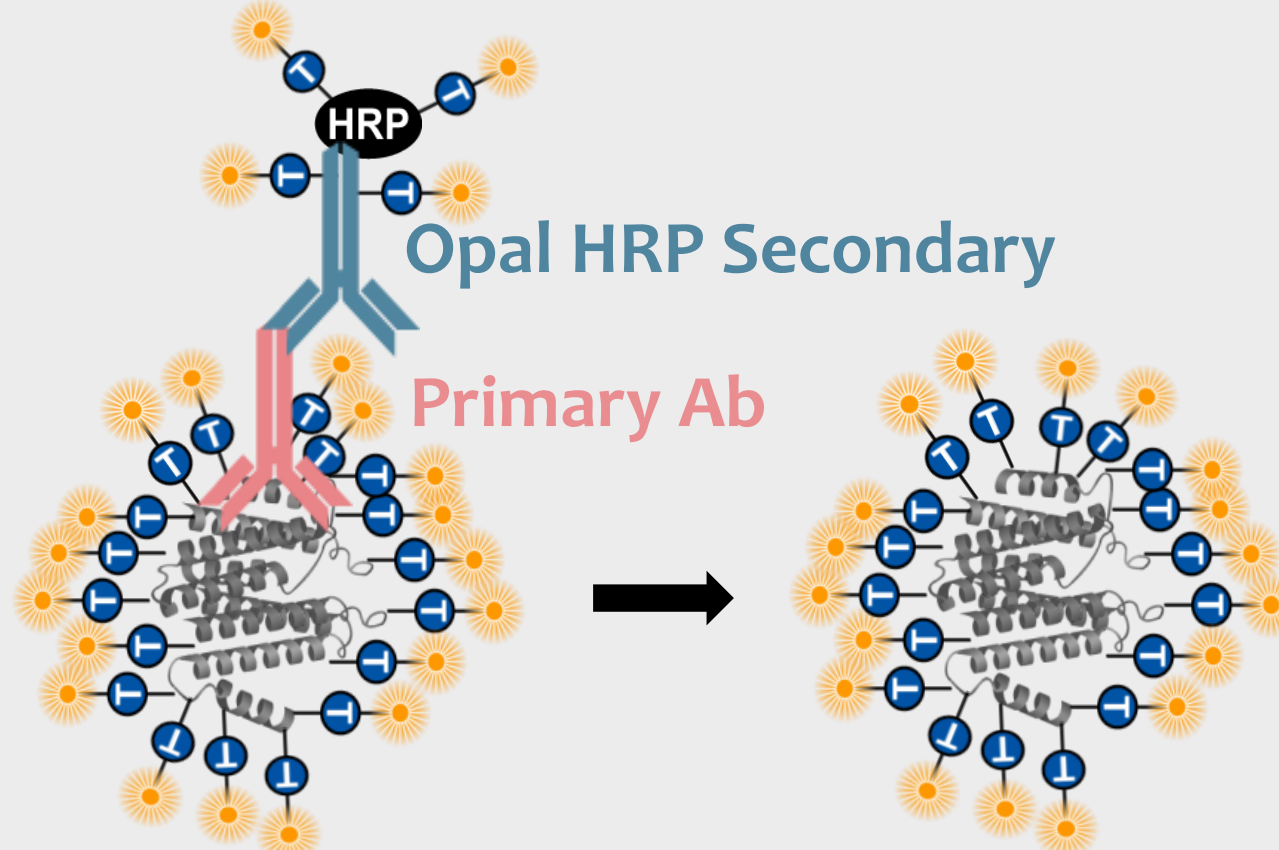


**Fig 1. Opal™ Detection.** The Opal HRP polymer amplifies IHC detection by covalently depositing multiple Opal fluorophores near the detected antigen. Then, antibodies are stripped to allow for sequential labeling of multiple markers.

### Multispectral Imaging and Unmixing

With Vectra Polaris®, different combinations of agile LED bands, bandpass excitation filters, bandpass emission filters, and a liquid crystal tunable filter (LCTF) are used to select narrow spectral bands for imaging. Images from up to 40 narrow spectral bands are layered into a single multispectral image.

**Fig 2. Unmixing in inForm® software.** (A) A multispectral image cube where each layer is taken at a different spectral range. (B) Signals from all spectral layers are linearly unmixed using pure emission spectra for each fluorophore. Intensity values in the resulting unmixed image (C) are directly related to the amount of each fluorophore present in each pixel.

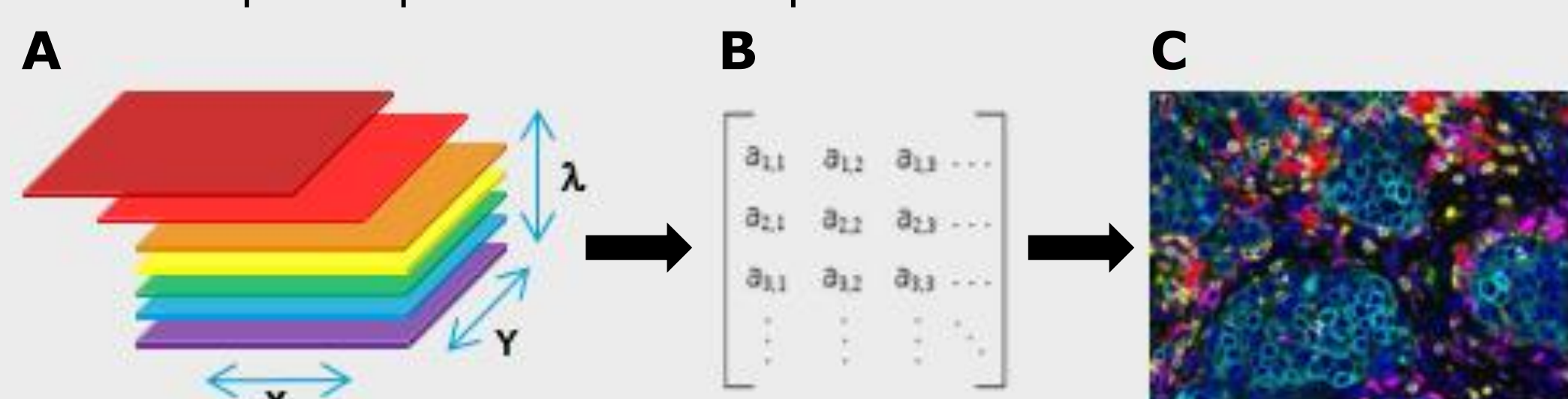


## 3 7-color Immuno-oncology Panel


FFPE tissue sections were stained with a 6-plex, 7-color Opal™ Polaris panel. The 6 markers in the panel combine to identify at least 12 different phenotypes of interest outlined below.

### Phenotype Combinations with 7-color Panel

Panel: PD-L1, PD-1, CD8, Tumor, DAPI, FoxP3, CD68

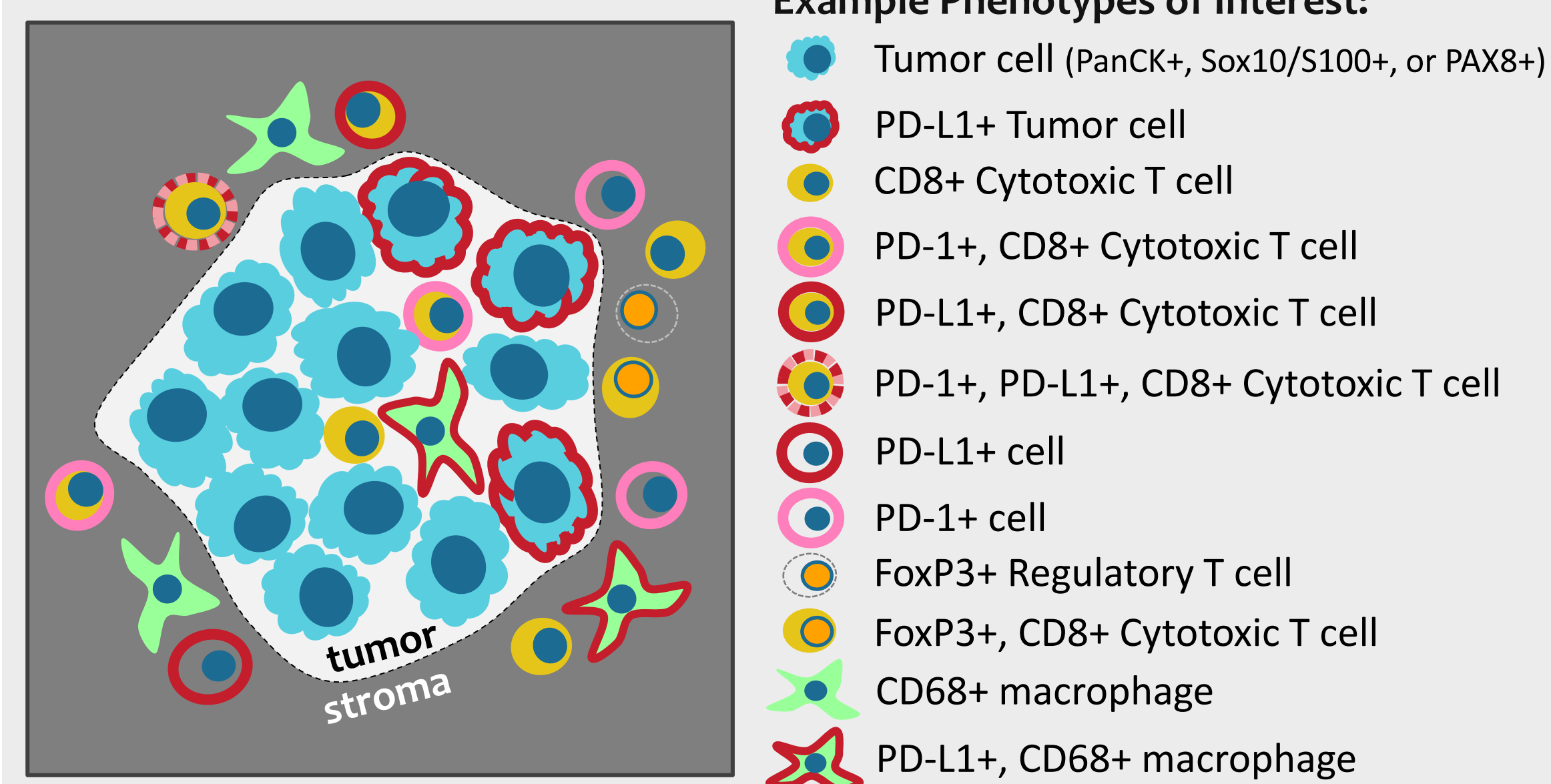


### Unmixed 6-plex, 7-color Multispectral Scan

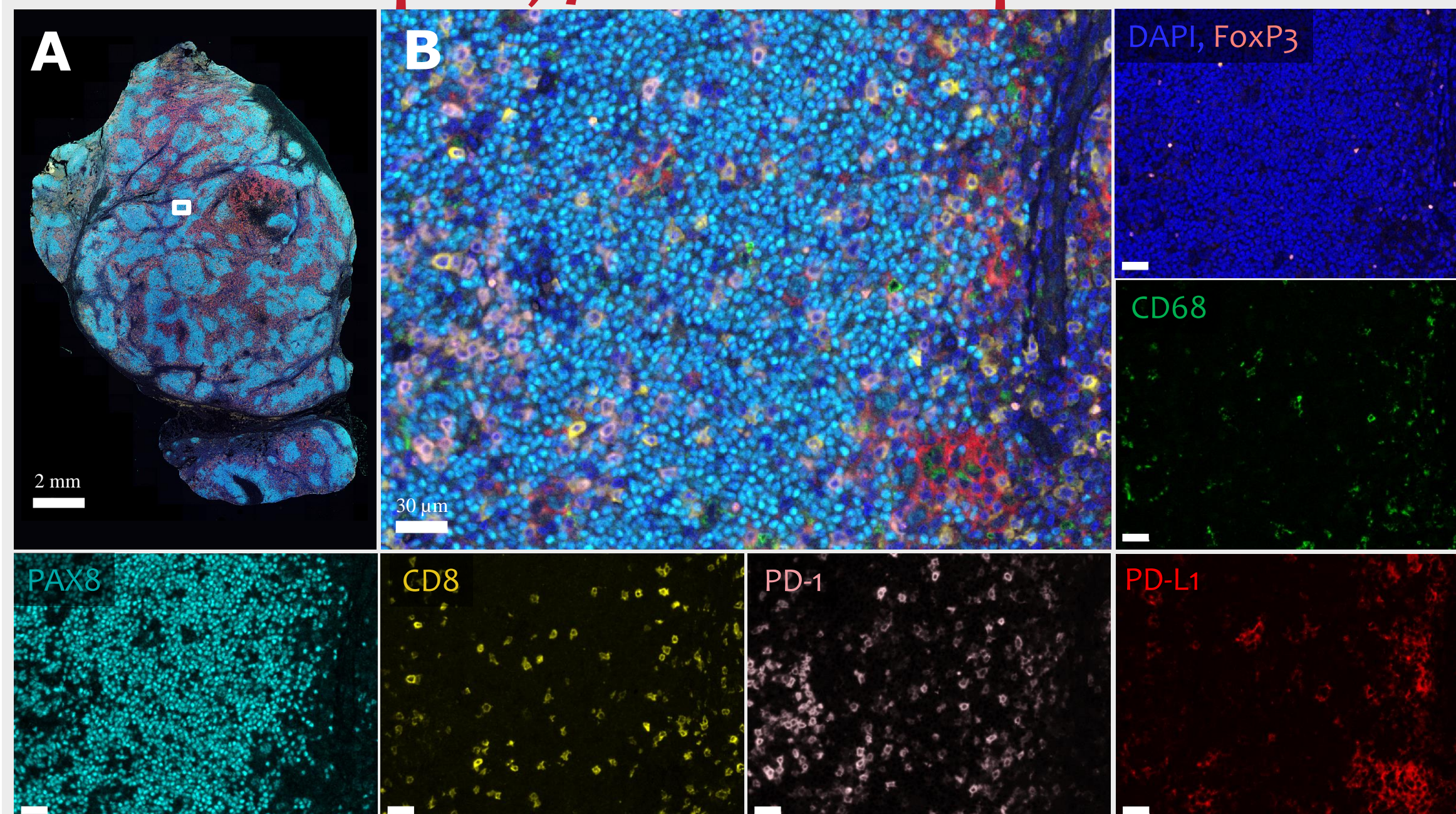

**Fig 3. Multispectral scan of human lymphoma tissue stained with 7-color Opal™ panel:** DAPI, PAX8 (Opal 690), CD8 (Opal Polaris 480), PD-1 (Opal Polaris 780), PD-L1 (Opal 520), CD68 (Opal 620), FoxP3 (Opal 570). (A) Whole slide unmixed view from Phenochart™. (B) View of region outlined by white rectangle in A. Remaining panels show separate planes that comprise B (scale bar, 30 μm).


## 4 Unmixing Improves Sensitivity

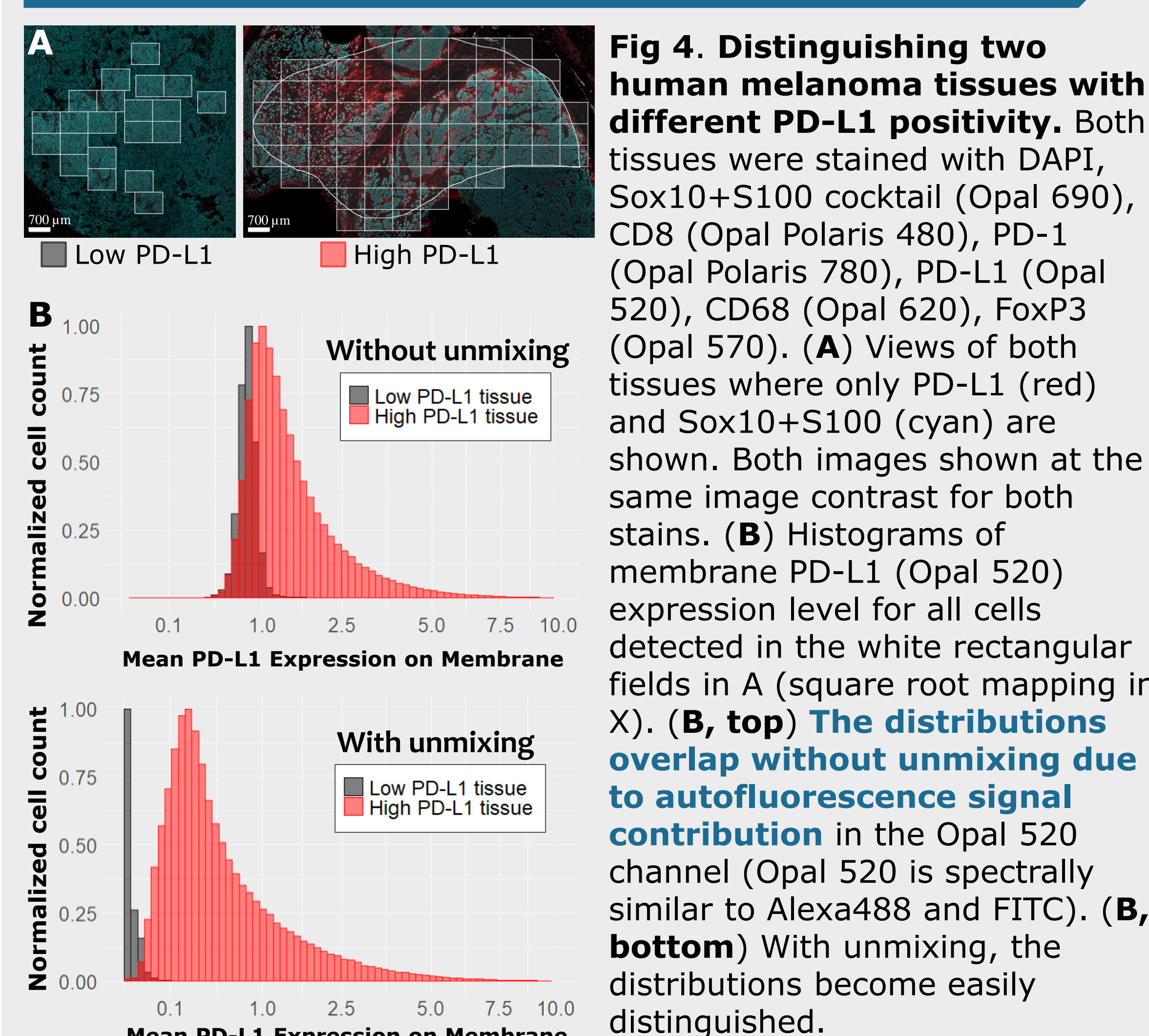

## 5 Characterizing the Tumor and Tumor Microenvironment


### Visualizing Heterogeneity within a Tissue Section

Unmixed images were processed with inForm® software to phenotype each marker in the panel. Resulting phenotype data can be plotted as heatmaps of cell density (cells/mm²). These density maps reveal heterogeneity across the section that is not immediately apparent from viewing the original image.

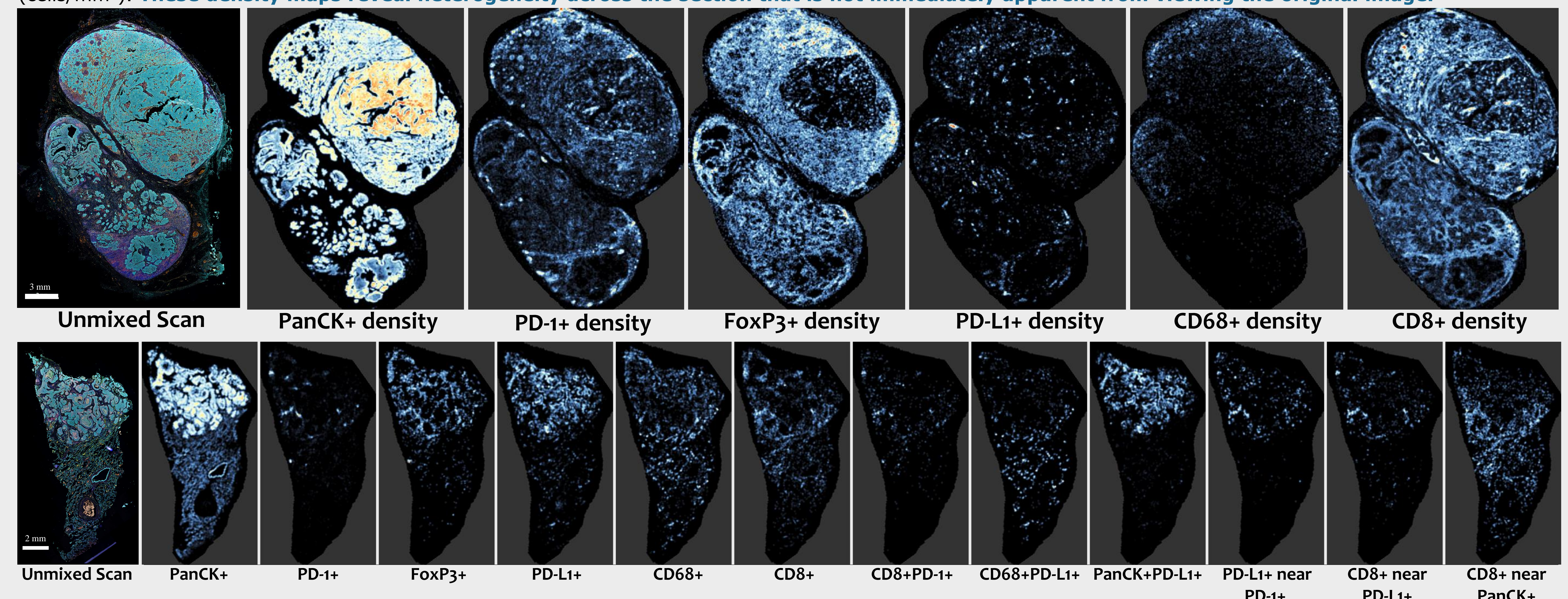


**Fig 5. Cell density maps of a head and neck cancer tissue (top) and lung cancer tissue (bottom).** 7-color Opal™ panel: DAPI, PanCK (Opal 690), CD8 (Opal Polaris 480), PD-1 (Opal Polaris 780), PD-L1 (Opal 520), CD68 (Opal 620), FoxP3 (Opal 570). Left-most panel is the unmixed multispectral scan. Remaining panels show heatmaps of cell density for different phenotypes (listed below each panel). Each panel has a different internal color scale. 'Near' indicates within 30 microns.

### Mapping Organization within the Tumor Microenvironment

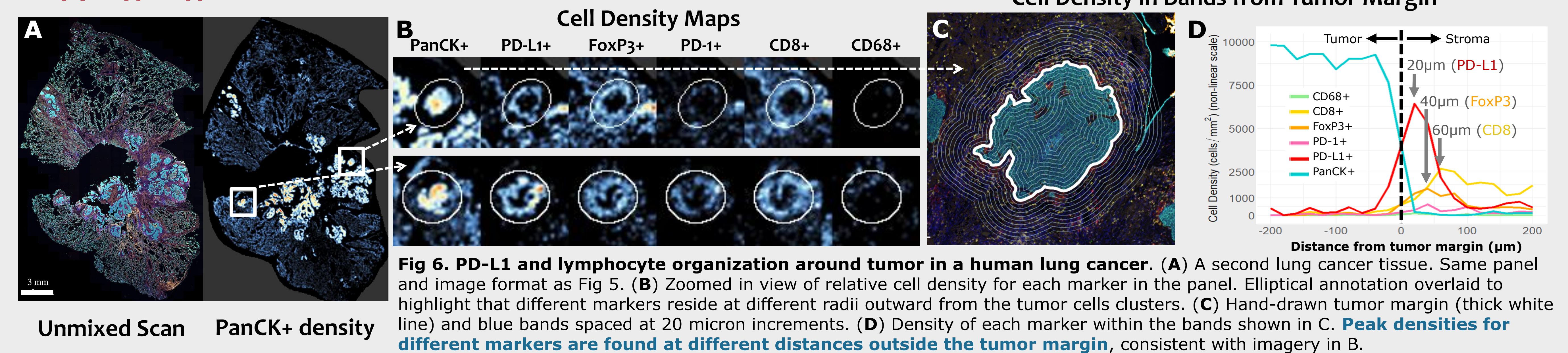

### Visualizing Heterogeneity Along a Tumor Margin

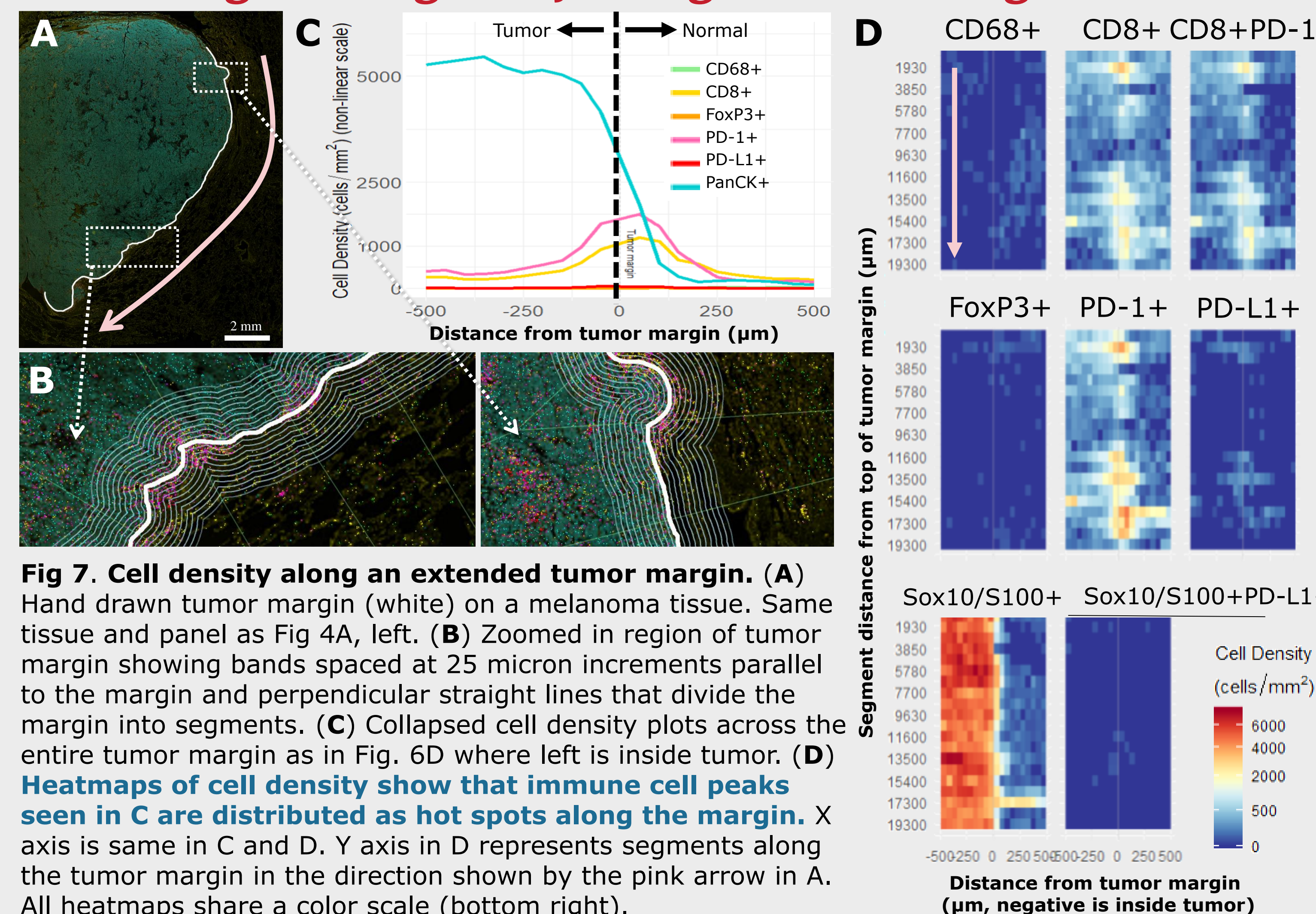

## 6 Conclusions

Spatial analyses of MOTIF™ multispectral scans showed wide variety among the tumor sections and within individual tumor sections. We observed differences in the overall density of individual markers, and further distinctions when looking at the local proximities of marker combinations that could not be observed with a conventional 3- or 4-color slide scan.

Furthermore, when compared with results from analogous scans analyzed without spectral unmixing, the MOTIF™ imagery showed a marked improvement in the dynamic range of detected positive cells.

These tools for spatial analyses of MOTIF™ high-throughput 7-color multispectral slide scans provide translational researchers the ability to probe a whole slide section with the necessary breadth and depth to describe the complex heterogeneity of the tumor microenvironment.

1. <https://akoyabio.github.io/phenoptReports/>  
2. RStudio Team (2015). RStudio: Integrated Development for R. RStudio, Inc., Boston, MA URL <http://www.rstudio.com/>.  
3. R Core Team (2019). R: A language and environment for statistical computing. R Foundation for Statistical Computing, Vienna, Austria. URL <https://www.R-project.org/>.  
4. Bankhead, P. et al. QuPath: Open source software for digital pathology image analysis. Sci. Rep. 7, 16878 (2017) <https://www.nature.com/articles/s41598-017-17204-5>